
## Instruction Manual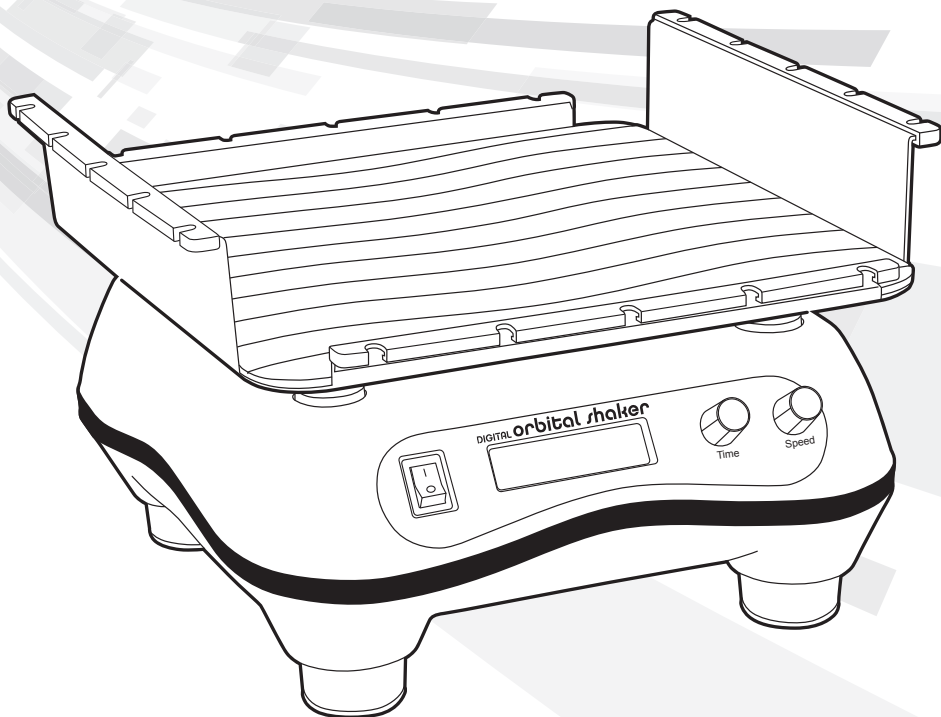

Instruction Manual.....	2
Manuel d'instructions.....	12
Manual de Instrucciones .....	22
Bedienungsanleitung .....	32

# Contents

Safety Precautions .....	2
Technical Data .....	3
Disposal Responsibilities .....	4

## Safety Precautions

This orbital shaker is a benchtop shaker designed for laboratories in the bioscience, physical science, and chemistry fields. It is NOT intended for use in CO<sub>2</sub> environments. Intended to mix solutions in various labware vessels. It may be involved with biohazardous and/or hazardous materials. This instruction manual can not address all safety hazards. It is the responsibility of the user to consult and observe all health and safety precautions and to assess the instrument's suitability to the task.

## Caution!

Not intended for CO<sub>2</sub> environments. Before starting the orbital shaker for the first time, please read this instruction manual carefully. This manual contains important information, safeguards, and operating instructions. Improper use of the instrument voids the warranty.

1. Before each use: inspect the unit for damage (broken or cracked housing, loose fasteners, damage from solvents, etc.) before connecting the power cord. Do not use the unit if any damage is found.
2. Never use this instrument in any manner inconsistent with these instructions.
3. Observe general laboratory safety precautions and regulations.
4. Do not modify construction and/or assembly of shaker.
5. Do not lift shaker by platform.
6. The platform must always be securely fastened.
7. All shaking vessels or containers must be secured to platform.
8. Do not use a damaged power adapter, power cord, platform, or bungee cords.
9. Use a properly grounded electrical outlet of correct voltage and current handling capacity.
10. Place the unit on a clean, sturdy, level surface at least 12" (30cm) away from the edge of the counter and other moving equipment. Make sure the AC power adapter and cords are clear of hot surfaces and other hazards. Ensure no person or hazardous substances are within a 12" (30cm) zone around the unit.
11. Stand clear of equipment when it is operating. Ensure that parts of the body, hair, or clothing can not be trapped by the motion of the unit.
12. Do not put hands in the platform area unless the platform is completely stopped.
13. Do not move unit while the shaker is shaking.
14. Do not immerse unit in water. Injury due to shock or fire may result.
15. The platform must be loaded symmetrically. Operating this unit with an unbalanced platform will cause severe vibration and damage, which is not covered by warranty.
16. Do not fill vessels or containers while attached to platform. Spilled liquid may harm unit.
17. Seal vessels or close containers before starting to shake. Open vessels or containers can spill contents during shaking and damage the unit.
18. Follow all precautions from the Material Safety Data Sheet (MSDS) for any reagent you use with this instrument.
19. Do not over extend bungee cord fasteners.
20. Use appropriate hand, eye, and face protection.

21. Do not use solvent or flammable liquids near this or other electrical equipment.
22. Do not shake flammable, explosive, or corrosive materials.
23. Do not operate in a hazardous or flammable environment.
24. Do not use any solvent on the unit that may attack plastic or cause damage.
25. Do not shake materials where the transfer of mechanical energy to glass could cause glass breakage.
26. If shaking action will result in the evolution of gases or fumes, carry out the operation in a well-ventilated laboratory hood.
27. Ensure the unit is protected from corrosion and mechanical damage. The unit must be cleaned with a pH-neutral cleaning liquid.
28. Always work in a manner, which endangers neither the user nor any other person.
29. Should the instrument fail to work to specification, immediately stop shaking. Clean and troubleshoot the instrument according to the instructions under "Trouble Shooting" before any further use of the instrument. Contact an authorized dealer or the manufacturer if trouble persists.
30. Disconnect from power supply before servicing or cleaning.
31. Do not stop platform while unit is turned ON. Doing so may cause the unit to fail and will void the warranty.
32. Only use original manufacturer's accessories (power adapter, platform, and bungee cords) and spare parts.
33. Repairs are to be performed by trained and authorized service personnel only. Contact Heathrow Scientific® LLC Repair Dept Technical Support.
34. Opening the instrument housing or improper use of the instrument voids the warranty. If there is a failure during the warranty period, contact Heathrow Scientific® LLC for warranty service.
35. Do not use units that have not been correctly installed or repaired.
36. Do not attempt to disassemble or modify.
37. Store at room temperature in a dry area. Do not expose to sunlight, moisture, or extreme temperatures or conditions for prolonged periods of time.

# Specifications

## Environmental Ratings

**Note:** For Indoor Use, Not for use in CO<sub>2</sub> environments

**Altitude:** up to 2,000m

**Temperature:** 2°C to 40°C

**Maximum Relative Humidity:** Non-condensing, 80% for temperatures up to 31°C decreasing linearly to 50% Relative Humidity at 40°C

### Storage Conditions:

Storage Temperature: 20°C - 55°C

Relative Humidity: <=90% noncondensing

## Technical Features

**Speed:** 20 to 300 rpm

**Orbit diameter:** 3/4" (19mm)

**Max load:** 4.5kg (10lbs)

**Timer:** 0 to 74 hours 59 minutes (4499 minutes) or continuous

### Dimensions (LxWxH):

Platform: Approx. 280mm x 280mm (11" x 11")

Device Overall: Approx. 315mm x 305mm x 192mm (12.38" x 12" x 7.56")

**Weight:** Approx. 9.80kg (21.6lb)

### Electrical:

#### Power Adapter Rating:

Input: 100-240VAC 50/60Hz 1.0A Comes with four interchangeable plugs for UK, Continental Europe/Israel, North America/Japan/Taiwan, Australia.

Output: 12VDC 2.2A

#### Shaker Rating:

Input: 12VDC

Current (under load) 1.5A

Power: 18W

## EC Declaration of Conformity

**Manufacturer:** Heathrow Scientific® LLC, 620 Lakeview Parkway, Vernon Hills, IL 60061 USA

**European Contact:** Emergo, Europe, Princessegracht 20, 2514AP, The Hague, The Netherlands

**Product:** Digital Orbital Shaker

**Product Model No(s):** 120460

This unit has been constructed and conforms to the following:

### EMC Standards and Directives:

EN61326-1:2013-b

FCC Part 15, Subpart: B: 2015

ICES-003, Issue 6

2014/35/EU (Low Voltage Directive) 2012/19/EU (WEEE)

RoHS3(EU 2015/863) and 2003/11/EC

### Safety Standards:

IEC/EN 61010-1:2010 ed. 3 + c1, c2

UL Std. 61010-1:2012 ed.3 + R:29 Apr2016

CAN/CSA C22.2 No. 61010-1-12:2012 ed.3+u1+u2(R2017)

Signed:

Gary Kamees

Title: Vice President Product Development/Manufacturing

Location: Vernon Hills, IL

Date of Issue: June 2022

## UKCA Declaration of Conformity

**Manufacturer:** Heathrow Scientific® LLC, 620 Lakeview Parkway, Vernon Hills, IL 60061 USA

**Product:** Digital Orbital Shaker

**Product Model No(s):** 120460

This unit has been constructed and conforms to the following:

### Directives and Regulations:

Electrical Equipment (Safety) Regulations 2016

The Restriction of the Use of Certain Hazardous Substances in Electrical and Electronic Equipment Regulations 2012

The Electromagnetic Compatibility Regulations 2016

### EMC Standards and Directives:

EN61326-1:2013-b

### Safety Standards:

UL Std. 61010-1:2012 ed.3 + R:29 Apr2016

CAN/CSA C22.2 No. 61010-1-12:2012 ed.3+u1+u2(R2017)

Signed:

Gary Kamees

Title: Vice President Product Development/Manufacturing

Location: Vernon Hills, IL

Date of Issue: October 2022

This mark is the confirmation that the unit conforms to the EU guidelines and has been tested according to the Electromagnetic Compatibility Directive 2014/30/EU and to the Low Voltage directive 2014/35/EU per the above EU Declaration of Conformity by an independent internationally recognized testing laboratory.

Please note that changes or modifications of this product is not expressly approved by the party responsible for compliance could void the user's authority to operate the equipment.

Note: this equipment has been tested and found to comply with the limits for a Class B digital device, pursuant to Part 15 of the FCC rules. These limits are designed to provide reasonable protection against harmful interference in a residential installation. This equipment generates, uses and can radiate radio frequency energy and, if not installed and used in accordance with the instructions, may cause harmful interference to radio communications. However, there is no guarantee that interference will not occur in a particular installation.


If this equipment does cause harmful interference to radio or television reception, which can be determined by turning by turning the equipment off and on, the user is encouraged to try to correct the interference by one or more of the following measures:

- Reorient or relocate the receiving antenna.
- Increase the separation between the equipment and receiver.
- Connect the equipment into an outlet on a circuit different from that to which the receiver is connected.
- Consult the dealer or an experienced radio/TV technician for help.

## QUICK START Guide

1. Read and understand all safety precautions and warnings first.
2. Make sure the specifications of this device are suitable for the application.
3. Switch power switch to OFF position.
4. Place the unit on a clean, sturdy, level surface at least 12" (30cm) away from the edge of the counter, table, or bench-top and other moving equipment. Ensure that no person or hazardous substances are within a 12" (30cm) zone around the units.
5. Connect the proper plug adapter to the AC power adapter, plug the power adapter into a wall outlet and the power cord into the rear of the device. Make sure the AC power adapter and cords are clear of hot surfaces and other hazards.
6. Secure vessel(s) or container(s) to be shaken to platform with included elastic cords.
7. Turn switch to ON position.
8. Rotate speed control to desired speed (displayed on screen). Press speed control knob to start unit.
9. Press speed control button to set timing or to stop shaking.
10. Adjust timer knob, press for continuous use.
11. To stop the device, press the speed control knob.

# Components

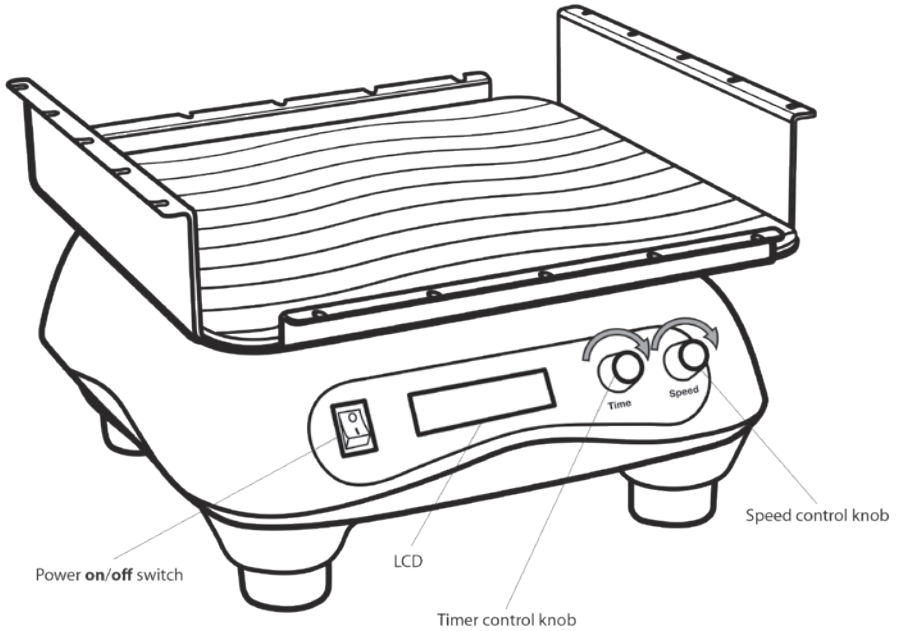## Disposal Responsibilities:

The equipment you purchased may contain hazardous substances that could impact the environment. Per regulations on electronic devices in the European Community, you must use the appropriate disposal systems to avoid exposure of these substances to the environment. The disposal systems will reuse or recycle hazardous materials from your equipment responsibly.

The crossed-out wheeled bin symbol invites you to use those systems.

If you need more information on the collection, reuse and recycling systems, please contact your local or regional waste administration. You can also contact us for more information on the environmental performance of our products.

# Getting Started

1. Remove the contents from the package, examining them carefully for breakage, defects, or missing parts. Save the packaging.

## **WARNING!**

**Do not lift shaker by the platform.**

2. Place shaker on a sturdy, level surface away 30 cm from the edge of the counter and other moving equipment. Allow 8 cm (3" ) of clearance around the unit for platform movement, accessory attachments, and user convenience. Make sure the power cord is clear of hot surfaces, sharp edges, and other hazards.

## **CAUTION!**

**If the top platform becomes jammed, it will not rotate. Keep platform free and clear of obstacles.**

3. Verify that the power switch is in the OFF position. Plug the power adapter cord into the rear of the shaker. Select the appropriate power plug attach to power adapter, and plug into an appropriate, grounded outlet.

## **WARNING!**

**Do not operate shaker with a damaged electrical cord.**

Note: — Leave shaker disconnected when not in use.

4. Turn power switch ON.
5. Rotate speed control knob counter clockwise until display reads 20 rpm. Press speed control knob once to turn shaker on. If there is a smooth whirring sound and the platform accelerates with little or no vibration, Digital Orbital Shaker is ready to use. If there are loud or unusual sounds, or excessive vibration, DO NOT OPERATE. Contact Heathrow Scientific® LLC Technical Support.
6. Before filling a container or vessel, it is recommended to test stability. Secure vessel on platform in middle position.
7. Press speed control knob once to turn on shaker. Make sure the vessel will not tip, teeter, walk, or slide to assure safe mixing.
8. Fill vessels for shaking. Vessels should contain a stopper or be covered to prevent spills.
9. Secure all vessels.

Notes: — Do not operate with an unbalanced load.

— Partial loads should be centered to evenly distribute weight over bearings for longer shaker life.

## To Set the Speed Function

1. Rotate the speed control knob clockwise to increase the speed and counterclockwise to decrease speed. The display will indicate the rpm set.
2. Press the speed control knob to start shaking or set timing function (see instructions below).
3. The speed control optical encoder provides instant readout of the actual platform rotation speed.

## To Set the Timer Function

1. Rotate the timer control knob clockwise to adjust the operating time. The display will indicate the minutes set.
2. Press the speed control knob to start shaking or change speed setting (see instructions above).
3. The speed control optical encoder provides instant readout of the actual platform rotation speed.
4. Unit will beep once programmed time has expired. Platform will stop shaking.
5. Remove vessels once platform has completely stopped.

## To Set the Continuous Timing Function

1. Press the time control knob to set for continuous use.
2. Press the speed control knob to start shaking or change speed setting (see instructions above).
3. The speed control optical encoder provides instant readout of the actual platform rotation speed.

## To Stop Shaking Function

1. Press the speed control knob to stop shaking.
2. Remove vessels once platform has completely stopped.

Note: — After an interruption in the power supply during shaking operation, the shaker will not start running again by itself.

# Maintenance

## **WARNING!**

**Disconnect plug from electrical outlet before attempting any maintenance.**

The shaker is normally maintenance-free. Clean unit only when it is not plugged into an electrical outlet. When necessary, the housing, platform, and mat can be wiped using a damp cloth and a mild, non-corrosive detergent. The mat must remain clean to maintain its friction-gripping quality.

## **WARNING!**

**The housing and platform must be cleaned with a neutral cleaning liquid.**

Note: — Ensure all parts are thoroughly dry prior to operation.

# Trouble Shooting

## **WARNING!**

**Disconnect plug from electrical outlet before attempting any maintenance.**

<b>Trouble</b>	<b>Possible Cause</b>	<b>Solution</b>
Screen says: *MOTOR OVERLOAD* PRESS RPM START	The top platform has jammed.	Clear obstacles. Press RPM knob.

If the above recommendation does not solve current problems, contact your authorized dealer or manufacturer directly.

# Return for Repair

## **IMPORTANT:**

**Transporting hazardous materials without a permit is a violation of U.S. federal law.**

Heathrow Scientific® LLC will not accept any shaker that is not appropriately cleaned and decontaminated. In the unlikely event of repair, or when damage to the unit necessitates return, contact Heathrow Scientific® LLC and obtain return authorization before sending your Digital Orbital Shaker for service.

# Warranty

## **LIMITED WARRANTY**

**Heathrow Scientific® LLC warrants that the shaker will be free from defects in workmanship and material for 5 years from the date of purchase.**

If you believe that there is a defect in the product, you must, during the warranty period, notify Heathrow Scientific® LLC, provide proof of purchase, and return the product to Heathrow Scientific® LLC with a Return Authorization form. To obtain a Return Authorization form, please call +1-847-816-5070.

If Heathrow Scientific® LLC is properly notified and, after inspection, confirms that there is a defect and the warranty period has not expired, Heathrow Scientific® LLC will repair, modify, or replace the product, at its sole option, at no charge.

OTHER THAN THIS LIMITED WARRANTY, HEATHROW SCIENTIFIC® LLC MAKES NO WARRANTIES, EXPRESS OR IMPLIED, WITH RESPECT TO THE QUALITY OR PERFORMANCE OF THE PRODUCT, INCLUDING BUT NOT LIMITED TO THE IMPLIED WARRANTIES OF MERCHANTABILITY AND FITNESS FOR PARTICULAR PURPOSE WHICH ARE HEREBY DISCLAIMED AND EXCLUDED. HEATHROW SCIENTIFIC® LLC WILL IN NO EVENT BE LIABLE FOR ANY LOSS OF USE, LOSS OF PROFITS, CONSEQUENTIAL, SPECIAL, EXEMPLARY OR PUNITIVE DAMAGES.

THIS WARRANTY DOES NOT COVER:

- ANY DEFECT OR DAMAGE CAUSED BY IMPROPER OR UNREASONABLE USE OF THE PRODUCT. (SHAKER IS DESIGNED FOR USE ONLY BY TRAINED LABORATORY TECHNICIANS. USE BY ANYONE ELSE WILL VOID THIS WARRANTY.)

- ANY PRODUCT THAT HAS BEEN, IN HEATHROW SCIENTIFIC® LLC'S SOLE JUDGEMENT, TAMPERED WITH, ALTERED, OR REPAIRED BY ANYONE OTHER THAN HEATHROW SCIENTIFIC® LLC.
- ANY PRODUCT THAT IS INOPERATIVE BECAUSE OF: (a) WEAR OCCASIONED BY USE, (b) NEGLIGENCE, (c) ACCIDENT, (d) INCORRECT MAINTENANCE, OR (e) USE UNDER ABNORMAL CONDITIONS OF TEMPERATURE, DIRT OR CORROSION, OR USE WITH ABRASIVE OR CORROSIVE MATERIALS.
- ACCESSORY PARTS, SUCH AS RUBBER PARTS, THAT ARE DAMAGED BY LIQUIDS OR MISUSE.

IN NO EVENT WILL HEATHROW SCIENTIFIC® LLC'S OBLIGATION UNDER THIS WARRANTY EXCEED THE PRICE OF THE PRODUCT.

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**HEATHROW SCIENTIFIC, LLC**