High-Speed Mini-Centrifuge

Instruction Manual	2
Manuel d'instructions.....	15
Manual de Instrucciones	29
Bedienungsanleitung.....	43

620 Lakeview Parkway • Vernon Hills, IL 60061
Phone: 1-847-816-5070 • Fax: 1-847-816-5072
Email: info@heathrowscientific.com
Web: www.heathrowscientific.com

Safety Precautions

Caution! Read entire manual and become familiar with the operation of this device before using it. If you have any questions, contact Heathrow Scientific LLC.

Your mini-centrifuge is designed for laboratories in the bioscience, medical and chemistry fields. This instruction manual cannot address all safety hazards. It does not have BIOSEALS and is not intended for use with radioactive materials. It is the responsibility of the user to consult and observe all health and safety precautions and to assess the instrument's suitability to the task.

CAUTION

Before starting the mini-centrifuge for the first time, please read this instruction manual carefully. This manual contains important information, safeguards, and operating instructions.

1. Never use this product in any manner inconsistent with these instructions.
2. This product is intended to be used indoors only.
3. This product is intended for separating aqueous solutions in approved test tubes compatible to the RCF of this unit.
4. The product is not intended to centrifuge very dense materials. Do not use materials with a density greater than 1.2 g/cm³ (1.2 g/mL).
5. Only use rotors supplied by Heathrow Scientific specifically made for this product. These rotors are rated for the maximum speed of the product, 12 500 rpm
6. Observe general laboratory safety precautions and regulations when using this product.
7. Do not attempt to operate with the cover removed or open.
8. The rotor and lid must always be securely fastened when in operation.
9. Always seat the rotor completely onto the shaft, then tighten the rotor nut.
10. Do not use damaged rotors.
11. Do not put hands onto the rotor unless the rotor is completely stopped.
12. Do not move unit while the rotor is spinning.
13. Do not immerse the product in water. Injury due to shock or fire may result.
14. The rotor must be loaded symmetrically. Operating this unit with an unbalanced rotor will cause severe vibration and damage, which is not covered by warranty.
15. Do not fill tubes while inserted in rotor. Spilled liquid may harm unit.
16. Close tube lids before starting centrifuge. Open lids can be torn off during centrifugation and damage the unit. Open lids can cause fluid to be dispersed inside the centrifuge.
17. Do not insert non-standard tubes or other foreign articles into the rotor.
18. Follow all precautions from the Material Safety Data Sheet (MSDS) for any reagent you use with this instrument.
19. Do not use solvent or flammable liquids near this or other electrical equipment.
20. Do not centrifuge materials that in combination or alone are flammable, explosive, corrosive or may cause any other hazard.
21. Do not operate in a hazardous or flammable environment.
22. Do not use any solvent on the unit that may attack plastic or cause cracks in the rotor.

23. Ensure the rotor is protected from corrosion and mechanical damage. The rotor must be cleaned with a pH-neutral cleaning liquid.
24. Always work in a manner which endangers neither the user nor any other person.
25. Should the instrument fail to work to specification, immediately stop centrifuging. Clean and troubleshoot the instrument according to the instructions under "Trouble Shooting" before any further use of the instrument. Contact an authorized dealer or the manufacturer if trouble persists.
26. Do not attempt to stop the rotor while unit is running. Doing so may cause the unit to fail and will void the warranty.
27. Only use original manufacturer's rotors, AC adapters, tube adapters and any other spare parts.
28. Repairs are to be performed by trained and authorized service personnel only. Contact Heathrow Scientific service.
29. Opening the instrument housing or improper use of the instrument voids the warranty. If there is a failure during the warranty period, contact Heathrow Scientific service for warranty service.
30. Do not attempt to use units that have not been correctly installed or repaired.
31. Do not attempt to disassemble or modify this product.
32. Store at room temperature in a dry area. Do not expose to sunlight, moisture, or extreme temperatures for prolonged periods of time.
33. Caution: Exposure to certain chemicals can damage product causing corrosion, crazing, and cracking of housing and components.
34. Before using any cleaning or decontamination methods except those recommended in this manual, please contact Heathrow Scientific to ensure that the proposed method will not damage the equipment.
35. If a single level of hazard protection is required, make sure to use containers that are tested to Risk Group II or above and that more than one level of protection is provided in the case of materials of a higher Risk Group.

WARNING

If the unit is not used as recommended by the manufacturer, the overall safety will be impaired.

WARNING

Check MSDS, wear required Personal Protective Equipment, and observe all applicable local and national regulations before dispensing and disposing of hazardous sample.

WARNING

The laboratory procedure should ensure that no person or hazardous substances are present within a 12" (30 cm) zone around the centrifuge when operating.

WARNING

This product does not contain bio-seals as per IEC/EN/CSA 61010-2-20 and cannot provide any level of containment in case of a spill or release of toxic, radioactive, or pathogenic micro-organisms thus these materials are not recommended to be used in this product.

Technical Data

Supplied Equipment

- 1 ea. Centrifuge
- 1 ea. Universal AC power adapter with 4 mains cables
- 1 ea. Standard Tube rotor (1.5 mL & 2.0 mL tubes) (12,500 rpm)
- 1 ea. Strip tube rotor (8 place 0.2 ml Strip Tubes or 32 individual 0.2 ml tubes) (12,500 rpm)
- 12 tube adapters for 0.2 mL tubes and 12 tube adapters for 0.5 mL tubes

Environmental Ratings:

Note: for indoor use

Altitude: up to 2 000 m

Temperature: 2° C to 40° C

Maximum Relative Humidity: 80% for temperatures up to 31° C decreasing linearly to 50% relative humidity at 40° C

Storage Conditions:

Storage temperature: 20° C – 55° C

Relative humidity: ≤ 90% noncondensing

Technical Features:

Capacity: (Max. liquid density: 1.2 g/cm³ or 1.2 g/mL):

Standard 12-position tube rotor: 24 mL Max.

Strip tube rotor: 6.4 mL Max,

Weight: Approx. 1.4 kg (3.1 lb) with 12-position rotor

Speed: 500 rpm - 12,500 rpm, Max RCF: 9 800 x g

Electrical:

Power Adapter rating:

Input: 100 – 240VAC, 50/60Hz, 2.0-1.0A

Output: 12VDC, 8.33A

Centrifuge:

Input: 12VDC,

Current (under load): Approx. 6.5A

Power: Approx. 78W

Positive Polarity

WARNING

ONLY use power cord and adapter supplied with unit!

EC Declaration of Conformity

Manufacturer: Heathrow Scientific® LLC, 620 Lakeview Parkway, Vernon Hills, IL 60061 USA

European Contact: Emergo, Europe, Princessegracht 20, 2514AP, The Hague, The Netherlands

Product: Gusto® High Speed Mini Centrifuge

Product Model No(s): HS10050, HS10061

This unit has been constructed and conforms to the following:

EMC Standards and Directives:

EN 61326-1:2013 (IEC61326-1:2012)

FCC Part 15, Subpart B

ICES-003, Issue 6

2014/35/EU (Low Voltage Directive)

2012/19/EU (WEEE)

RoHS3 (EU 2015/863)

Safety Standards:

EN/IEC61010-1ed.3.1, 2017

EN/IEC61010-2-020 ed.3, 2016

UL Std. 61010-1 ed.3.1, 2017

UL Std. 61010-2-020 ed.3, 2016

CAN / CSA C22.2 No. 61010-1-12 (R2017)

CAN/CSA C22.2 No. 61010-2-020 (R2017)

Signed:

Gary Kamees

Title: Vice President Product Development/Manufacturing

Location: Vernon Hills, IL

Date of Issue: June 2022

UKCA Declaration of Conformity

Manufacturer: Heathrow Scientific® LLC, 620 Lakeview Parkway, Vernon Hills, IL 60061 USA

Product: Gusto® High Speed Mini Centrifuge

Product Model No(s): HS10050, HS10061

This unit has been constructed and conforms to the following:

Directives and Regulations:

Electrical Equipment (Safety) Regulations 2016

The Restriction of the Use of Certain Hazardous Substances in Electrical and Electronic Equipment Regulations 2012

The Electromagnetic Compatibility Regulations 2016

EMC Standards:

EN 61326-1:2013 (IEC61326-1:2012)

ICES-003, Issue 6

Safety Standards:

EN/IEC61010-1ed.3.1, 2017

EN/IEC61010-2-020 ed.3, 2016

UL Std. 61010-1 ed.3.1, 2017

UL Std. 61010-2-020 ed.3, 2016

CAN / CSA C22.2 No. 61010-1-12 (R2017)

CAN/CSA C22.2 No. 61010-2-020 (R2017)

Signed:

Gary Kamees

Title: Vice President Product Development/Manufacturing

Location: Vernon Hills, IL

Date of Issue: October 2022

NOTICE

Changes or modifications to the product not expressly approved by the manufacturer could void the user's authority to operate the equipment.

NOTICE

This product has been tested and found to comply with the limits of a Class A digital device pursuant to Part 15 of the FCC rules. These limits are designed to provide reasonable protection against harmful interference when the product is operating in a commercial environment. The product generates, uses, and can radiate radio frequency energy and if not installed and used in accordance with the user's manual may cause harmful interference to radio communications. Operation of the product in a residential area is likely to cause interference in which case the user will be required to correct the interference.

Disposal Responsibilities


The equipment you purchased may contain hazardous substances that could impact the environment. Per regulations on electronic devices in the European Community, you must use the appropriate disposal systems to avoid exposure of these substances to the environment. The disposal systems will reuse or recycle hazardous materials from your equipment responsibly.

The crossed-out wheeled bin symbol invites you to use those systems.


If you need more information on the collection, reuse and recycling systems, please contact your local or regional waste administration. You can also contact Heathrow Scientific for more information on the environmental performance of our products.

Display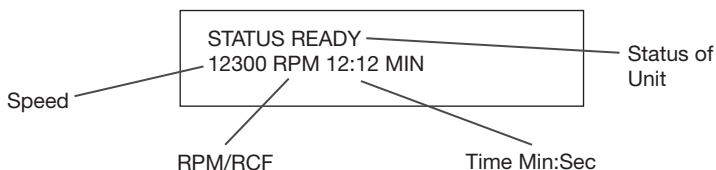


DISPLAY STATUS	
LID OPEN	- Lid is open
READY	- Unit ready to centrifuge
SPIN UP	- Cycle is starting and is increasing speed
SPINNING	- Running in Standard mode
SPIN DOWN	- Cycle is slowing down
STOPPED	- Cycle has stopped
COMPLETED	- Centrifuge cycle is finished
QUICK SPIN (flashes)	- Running in Quick Spin mode
ERROR	- An error has occurred (See error section)

FRONT PANEL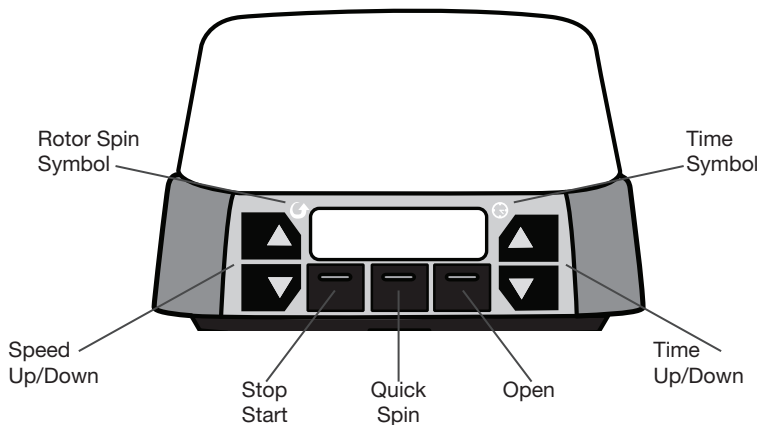

REAR PANEL

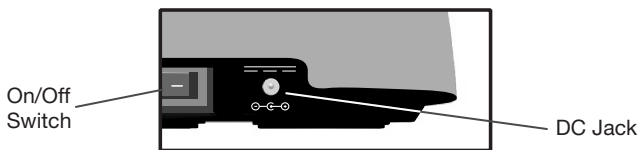Operation

- Remove the contents from the package examining them carefully for breakage, defects or missing parts.
- Place the product on a sturdy, level surface at least 300mm away from the edge of the counter and other moving equipment. Make sure the AC power adapter and cords are clear of hot surfaces and other hazards.
- Make sure the product has ventilation and is not encased in any material that will limit airflow. An overheating situation could occur.
- Verify that the power switch in the back of the unit is in the OFF position. Plug the DC cord into the rear of the unit and the AC cord into an approved outlet.
- **ONLY** use Heathrow Scientific supplied power adapter and cords.

Prior to using

- Before every use, inspect unit for cracks, missing pieces, and or any damage that would result in an impaired performance and notify manufacturer BEFORE use.
- Turn power switch on the rear ON. The unit will initialize.
- Press the OPEN button. The lid should pop up slightly. Lift the front of the lid gently up and backward until approximately 90 degrees vertical.
- Remove any remaining packaging materials.
- Verify that the rotor is installed correctly. (see section Rotor Installation and Removal) Do not run without the rotor installed.
- Close the lid.
- Press the START/STOP button (rotor should spin up to the preset speed). If there is a smooth whirring sound and the unit accelerates with little or no vibration, the mini-centrifuge is ready to use. If there are loud or unusual sounds, or excessive vibration, DO NOT OPERATE. Contact Heathrow Scientific service.
- At the set time, the rotor will spin down, stop and the unit will emit a beep when the cycle is complete.

General Operation

1. Turn power switch on the back of the unit ON. The unit will initialize.
2. To open lid, press the OPEN button. The lid should pop up slightly. Lift the front of the lid gently up and backward until approximately 90 degrees vertical.
3. On benchtop, prepare samples tubes in a rack so that tubes are filled to equal levels. Close tube lids.
4. Install tubes in the rotor in a balanced manner (see section - Balancing the Rotor).
5. Set the appropriate time by pressing the TIME UP/DOWN button.
6. Set the appropriate rpm by pressing the SPEED UP/DOWN button.
7. Place the standard rotor cover on the rotor to limit noise and aspiration of liquids that may escape from tubes.
8. Close Lid and press down until locked (Display status should read "READY").
9. Press START/STOP (rotor should spin up to the user set speed).
10. At a determined time the rotor will spin down and stop at the preset time.
11. The unit will then emit a beep when the cycle is complete.
12. To open lid, press the OPEN button.

Quick Spin Operation

1. Prepare the unit for centrifugation as in the General Operation (section above Steps 1-8).
2. Press and HOLD QUICK SPIN button (rotor should spin up to the set speed).
3. The timer will begin counting.
4. Once enough time has elapsed for your needs, release the QUICK SPIN button. The rotor will spin down, stop and the unit will emit a beep when the cycle is complete.
5. To open lid, press the OPEN button.

Other

Change Between RPM and RCF


- Prior to a spin cycle, press both the SPEED UP & DOWN buttons simultaneously.
- The speed setting will then convert to the other.

Quick Stop

- If for any reason you need to stop the cycle quickly, press and hold the START/STOP button. The cycle will stop more rapidly than normal. Caution The tube contents may be disturbed during the quick stop process.

Error

- If an error happens, the unit will beep and the display will indicate the error.
- Press Start/Stop to clear the error from the display.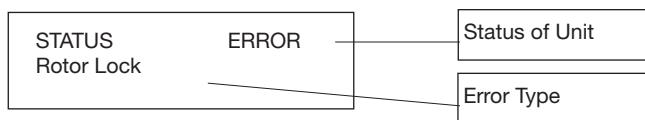

Error Status	Resolution
Motor Overload	<ul style="list-style-type: none">• Something is interfering with the rotor. Clear the rotor and reset.
User Stop	<ul style="list-style-type: none">• The operator has held down the START/STOP and implemented a quick stop.
Balance	<ul style="list-style-type: none">• Inspect the tubes for equal tube fill or improper placement. Once corrected, rerun.• If the balance error continues to happen, remove the tubes and determine if the balance error still persists with an empty rotor.• Inspect the rotor for improper installation.
Temperature	<ul style="list-style-type: none">• The unit has exceeded the normal operating temperature.• Turn off the unit and allow to cool.
Excessive Tilt	<ul style="list-style-type: none">• The unit has experienced a non-normal tilt event.• Make sure the unit is placed on a level surface. Once corrected, rerun
Lid Fail	<ul style="list-style-type: none">• The lid has opened during the cycle. Check for proper operation of the lid lock mechanism. The lid should stay locked during the entire cycle.
Rotor Lock	<ul style="list-style-type: none">• The unit has experienced a problem with the rotor.• Correct the rotor interference. Once corrected, rerun.

Trouble Shooting

Trouble	Resolution
No power present	<ul style="list-style-type: none">• Verify that the AC adapter is fully plugged into the wall and rear of unit.• Verify that the power switch is turned on.
Unit or display is not operating normally	<ul style="list-style-type: none">• Turn off the unit, wait 2 minutes, and turn power on.
Excessive vibration Excessive noise	<ul style="list-style-type: none">• Inspect the tubes for equal fill or improper placement.• Inspect the rotor for improper installation.• Remove the tubes and determine if the noise persists with an empty rotor.
Lid will not close	<ul style="list-style-type: none">• Verify that nothing is blocking the lid from fully closing.• Verify that nothing has fallen into the lock mechanism opening.
Lid will not open	<p>If for any reason you need to manually open the lid to access tubes due to an error or power loss, please perform the below:</p> <ul style="list-style-type: none">• Turn off the unit, remove power cord.• Make sure the rotor has stopped completely.• Use a thin rod and insert it into the opening on the bottom.• Press gently but firmly. You will feel a mechanical movement within the unit and the lid will release.• Remove the rod, set the unit on the feet.• Remove your tubes and reclose the lid.

WARNING


Spin balanced loads only!

Do not place the tubes asymmetrically nor load unequal volume tubes. Improper placement of tubes will lead to insufficient centrifugation and may cause serious injury or an accident.


BALANCING THE ROTOR

Tubes of equal weight and size should be placed opposite each other. Use additional sample or water in other tubes to provide a balanced rotor. Examples of proper sample balancing are illustrated next.

1.5 mL / 2.0 mL Circular Rotor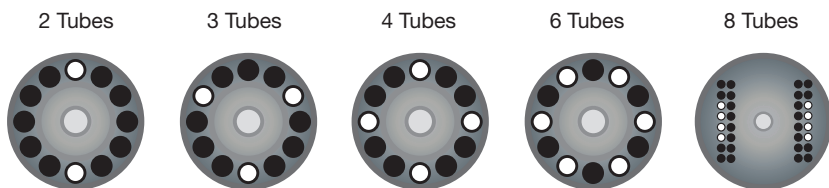

0.2 mL Tube Rotor

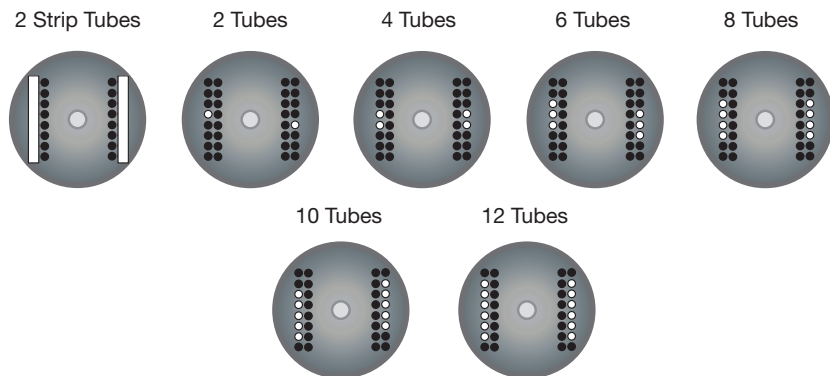WARNING

Do not attempt to open lid until rotor has completely stopped!
Avoid severe personal injury or property damage from moving parts.
Only use rotor compatible with the centrifuge.

Rotor installation and removal

To remove rotor:

- With one hand, grasp the rotor. With the other hand, unscrew center knob counter clockwise until it disengages from the motor shaft.
- Pull up on the center knob firmly. The rotor should easily lift off of the motor drive.

To install the rotor:

- Check to make sure the bands on the rotor are in the slots.
- Set rotor onto the motor drive and push down until seated.
- With one hand, grasp the rotor. With the other hand, screw the center knob clockwise until it engages the motor shaft. Keep turning the knob until it tightens.
- Do not overtighten the knob.

WARNING

Do not use damaged or cracked rotors!

Maintenance

Your mini-centrifuge is maintenance-free. Clean unit only when it is not plugged into an electrical outlet. When necessary, the housing and rotor can be wiped using a damp cloth and a mild, non-corrosive detergent.

WARNING

Do not use any solvent on the unit that may attack plastic or cause cracks in the rotor. Ensure the rotor is protected from corrosion and mechanical damage. The rotor must be cleaned with a neutral cleaning liquid.

If hazardous sample is spilled on the unit and or leaked into the unit, do not run the centrifuge unless it is safe to do so.

Dissassembly of Rotors for Cleaning

1. Remove rotor from unit
2. Remove screws from the bottom of the rotor
3. Clean as needed
4. Reassemble rotor, making sure that the alignment tab is engaged into the slot.
5. Test run reassembled empty rotor to ensure it is still balanced

NOTICE

Ensure all parts are thoroughly dried prior to operation.

Hazardous Chemicals:

1. Cleaning and decontamination may be necessary as a safeguard before laboratory centrifuges, rotors, and any accessories are maintained, repaired, or transferred. The user must follow any and all instructions provided in the manual pertaining to this process.
2. In the event hazardous material is spilt, the user is responsible for carrying out the appropriate decontamination procedure.
3. Before using any cleaning or decontamination methods except those recommended by the manufacturer, users should check with the manufacturer that the proposed methods will not damage the equipment.

Ordering Information

Mini-centrifuge with AC adapter, 2 rotors, rotor cover, and tube adapters.....	HS10050
Standard Rotor.....	HS100503
Strip Tube Rotor.....	HS100504
Rotor Knob.....	HS100505
Standard Rotor Cover.....	HS100506
Cord Set.....	HEA100502
Power Adapter.....	HEA120527
Tube Adapters (0.2 mL x6, 0.5 mL x6).....	x2 HEA120015

Return for Repair

IMPORTANT

Transporting hazardous materials without a permit is a violation of federal law.

Heathrow Scientific will not accept any product return that is not appropriately cleaned and decontaminated. In the unlikely event of repair, or when damage to the unit necessitates return, contact Heathrow Scientific, obtain return authorization, and complete a Heathrow decontamination form BEFORE sending your product for service.

Warranty

IN NO EVENT WILL HEATHROW SCIENTIFIC® LLC'S OBLIGATION UNDER THIS WARRANTY EXCEED THE PRICE OF THE PRODUCT.

Limited Warranty

Heathrow Scientific warrants that your mini-centrifuge will be free from defects in workmanship and material for 5 years from the date of purchase.

If you believe that there is a defect in the product, you must, during the warranty period, notify Heathrow Scientific, provide proof of purchase, and return the product to Heathrow Scientific with a Return Authorization form.

If Heathrow Scientific is properly notified and, after inspection, confirms that there is a defect and the warranty period has not expired, Heathrow Scientific will repair, modify, or replace the product, at its sole option, at no charge.

OTHER THAN THIS LIMITED WARRANTY, HEATHROW SCIENTIFIC MAKES NO WARRANTIES, EXPRESS OR IMPLIED, WITH RESPECT TO THE QUALITY OR PERFORMANCE OF THE PRODUCT, INCLUDING BUT NOT LIMITED TO THE IMPLIED WARRANTIES OF MERCHANTABILITY AND FITNESS FOR PARTICULAR PURPOSE WHICH ARE HEREBY DISCLAIMED AND EXCLUDED. HEATHROW SCIENTIFIC WILL IN NO EVENT BE LIABLE FOR ANY LOSS OF USE, LOSS OF PROFITS, CONSEQUENTIAL, SPECIAL, EXEMPLARY OR PUNITIVE DAMAGES.

THIS WARRANTY DOES NOT COVER:

- ANY DEFECT OR DAMAGE CAUSED BY IMPROPER OR UNREASONABLE USE OF THE PRODUCT. (THE MINI-CENTRIFUGE IS DESIGNED FOR USE ONLY BY TRAINED LABORATORY TECHNICIANS. USE BY ANYONE ELSE WILL VOID THIS WARRANTY.)

- ANY PRODUCT THAT HAS BEEN, IN HEATHROW SCIENTIFIC SOLE JUDGEMENT, TAMPERED WITH, ALTERED, OR REPAIRED BY ANYONE OTHER THAN HEATHROW SCIENTIFIC.
- ANY PRODUCT THAT IS INOPERATIVE BECAUSE OF: (a) WEAR OCCASIONED BY USE, (b) NEGLIGENCE, (c) ACCIDENT, (d) INCORRECT MAINTENANCE, OR (e) USE UNDER ABNORMAL CONDITIONS OF TEMPERATURE, DIRT OR CORROSION, OR USE WITH ABRASIVE OR CORROSIVE MATERIALS.
- ACCESSORY PARTS, SUCH AS RUBBER AND PLASTIC PARTS THAT ARE DAMAGED BY LIQUIDS OR MISUSE.

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All rights reserved. Printed in China.
©Heathrow Scientific® LLC