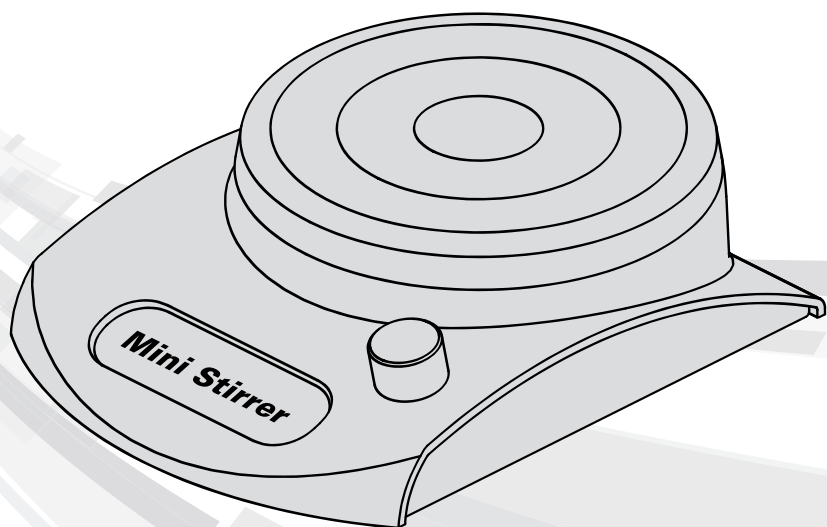
Instruction Manual

Instruction Manual	2
Manuel d'instructions	10
Manual de Instrucciones	18
Bedienungsanleitung	26

Safety Precautions

The "Mini Stirrer" is a small Magnetic Stirrer designed for laboratories in the bioscience, medical, and chemistry fields. This instruction manual cannot address all safety hazards. It is the responsibility of the user to consult and observe all health and safety precautions and to assess the instruments suitability to the task.

Caution!

Before operating the mini Stirrer please read this instruction manual carefully. This manual contains important information, safeguards, and operating instructions.

1. Always inspect the unit for damage (broken or cracked housing and base, loose fasteners, damage from solvents, etc.) before connecting the power cord and before each use. Do not use the unit if any damage is found.
2. Never use this product in any manner inconsistent with these instructions.
3. This product is intended to be used indoors only.
4. This product is intended for blending aqueous solutions in suitable approved containers.
5. Product is not suitable for use with high viscosity materials, some examples are; heavy weight oils, honey, or glycerin.
6. Use only the power adapter delivered with the product, use of other power adapters will void the warranty.
7. Carefully note position of container, efficiency is maximized when container and base are concentric, see molded in concentric circles on top of unit for various container sizes.
8. Observe general laboratory safety precautions and regulations.
9. Do not move the unit while it is operating.
10. Do not immerse the unit in water. Injury due to shock or fire may result.
11. Follow all precautions from the material safety data sheet (MSDS) for any reagents you use with this instrument.
12. Do not mix materials that are explosive.
13. Do not operate in a hazardous or flammable environment.
14. Do not use solvent or flammable liquids near this or other electrical equipment.
15. Do not use any caustic solvents on the unit that may damage the plastic housing.
16. Always work in a safe manner.
17. Opening the instrument housing or improper use of the instrument voids the warranty. If there is a failure during the warranty period, contact Heathrow Scientific LLC for warranty replacement.

18. Do not attempt to disassemble or modify this product.
19. Store at room temperature in a dry area. Do not expose to sunlight, moisture, or extreme temperatures for prolonged periods of time.

Warning!

If the unit is not used as recommended by the manufacturer, injury may result.

Warning!

Check MSDS, wear required Personal Protective Equipment, and observe all applicable local and national regulations before dispensing and disposing of hazardous samples.

Warning!

Do not use any solvent on the unit that may attack plastic. If hazardous sample material is spilled on the unit and or leaked into the unit, do not run the unit unless it is safe to do so.

Warning!

The laboratory procedure should ensure that no person or hazardous substances are present within a 12" (30cm) zone around the unit before operating.

Specifications

Environmental Ratings:

Note: for Indoor use

Altitude: up to 2 000m

Temperature: 2°C to 40°C

Maximum Relative Humidity: Non-condensing, 80% for temperatures up to 31°C decreasing linearly to 50% relative humidity at 40°C

Storage Conditions:

Storage Temperature: 20°C - 55°C

Relative humidity: ≤ 90% noncondensing

Technical Features:

Capacity:

Max Liquid volume: 1000 ml (1 Liter)

Dimensions:

180mm x 137mm x 50mm (LxWxH)

7.1" x 5.4" x 2.0" (LxWxH)

Weight:

Approx. 1 lb (.45 kg) not including power supply

Maximum Speed:

Unloaded Approx: 2400 RPM

Loaded Approx: 2000 RPM

Minimum Speed:

Unloaded Approx: 200 RPM

Speed Indication and Control Knob:

Located on front of unit

Electrical:

Power Adapter Rating

Input: 100-240VAC

50/60Hz

0.3A

Output: 12VDC 0.5A

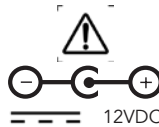
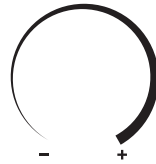
Device comes with interchangeable plugs for the UK, Europe, North America, and Australia.

Mini Stirrer Rating

Input: 12VDC

Current (under load): Approx. 0.4A

Power: Approx. 4.8W

This mark is the confirmation that the unit conforms to the EU guidelines and has been tested according to the following EU Declaration of Conformity.

Conforms to:
IEC/EN 61010-1:2010 ed.3 + c1, c2
ANSI/UL 61010-1:2012 ed.3 (R2016)
IEC 61010-2-051: ed.3
ANSI/UL 61010-2-051:2015 ed.3
CAN/CSA C22.2 No. 61010-1-12 (R 2017)
CAN/CSA C22.2 No. 61010-2-051:2015 ed.3

EMC Standards:
IEC 61326-1:2012 (EN 61326 1:2013)
FCC Part 15, Subpart B
ICES-003, Issue 6
2014/35/EU (Low Voltage Directive)
2012/19/EU (WEEE)
RoHS3 (EU 2015/863)

Intertek

EC Declaration Of Conformity

Manufacturer: Heathrow Scientific® LLC, 620 Lakeview Parkway, Vernon Hills, IL 60061 USA

European Contact: Emergo, Europe, Princessegracht 20, 2514AP, The Hague, The Netherlands

Product: Mini Stirrer

Product Model No(s): 120155, 120595

This unit has been constructed and conforms to the following:

EMC Standards and Directives:

IEC 61326-1:2012 (EN 61326 1:2013)

FCC Part 15, Subpart B

ICES-003, Issue 6

2014/35/EU (Low Voltage Directive)

2012/19/EU (WEEE)

RoHS3 (EU 2015/863)

Safety Standards:

IEC/EN 61010-1:2010 ed.3 + c1, c2

ANSI/UL 61010-1:2012 ed.3 (R2016)

IEC 61010-2-051: ed.3

ANSI/UL 61010-2-051:2015 ed.3

CAN/CSA C22.2 No. 61010-1-12 (R 2017)

CAN/CSA C22.2 No. 61010-2-051:2015 ed.3

Signed:

Gary Kamees

Title: Vice President Product Development/Manufacturing

Location: Vernon Hills, IL

Date of Issue: June 2022

UKCA Declaration Of Conformity

Manufacturer: Heathrow Scientific® LLC, 620 Lakeview Parkway, Vernon Hills, IL 60061 USA

Product: Mini Stirrer

Product Model No(s): 120155, 120595

This unit has been constructed and conforms to the following:

Directives and Regulations:

Electrical Equipment (Safety) Regulations 2016

The Restriction of the Use of Certain Hazardous Substances in Electrical and Electronic Equipment Regulations 2012

The Electromagnetic Compatibility Regulations 2016

EMC Standards:

IEC 61326-1:2012 (EN 61326 1:2013)

ICES-003, Issue 6

Safety Standards:

IEC/EN 61010-1:2010 ed.3 + c1, c2

IEC 61010-2-051: ed.3

CAN/CSA C22.2 No. 61010-1-12 (R 2017)

ANSI/UL 61010-1:2012 ed.3 (R2016)

ANSI/UL 61010-2-051:2015 ed.3

CAN/CSA C22.2 No. 61010-2-051:2015 ed.3

Signed:

Gary Kamees

Title: Vice President Product Development/Manufacturing

Location: Vernon Hills, IL

Date of Issue: October 2022

Operation:

1. Remove the contents from the package, examining them carefully for breakage, defects, or missing parts. Save the packaging.
2. Place Mini Stirrer on a sturdy, stable, level surface near a grounded electrical outlet away from the edge of the counter and other moving equipment. The surface should be clean. Allow sufficient clearance on all sides of the unit for proper ventilation and usage. Make sure the power cord is clear of hot surfaces and other hazards.
3. Place stir bar into bottom of empty container (see warning below).
4. Carefully fill container with all fluids to be mixed, note the level of the liquid which will rise to accommodate the spinning vortex of fluid when unit is powered. Be sure to leave sufficient room at the container top for this rise in fluid level.
5. Center the mixing container on top of the stirrer, noting and verifying the magnetic coupling of the stir-bar.
6. Turn control knob clockwise to turn on unit and continue steadily turning until desired mixing speed is obtained.
7. Turn off unit before removing mixing container.

Warning!

Do not drop stir bars into empty glass containers as the container may break.

Warning!

Do not plug other power supplies into stirrer unit, this will void warranty.

Warning!

Do not use if cord or power supply is damaged.

Maintenance

The Mini Stirrer is a maintenance free product. Clean unit only when it not plugged into an electrical outlet. When necessary, the complete outside of the unit can be wiped down using a damp cloth and a mild, non-corrosive detergent. Do not immerse the unit or the power adapter.

Trouble Shooting

Trouble	Possible Cause	Solution
Stir bar is not rotating	— Not properly plugged in	— Verify unit is plugged in
	— Unit turned off	— Turn unit on
	— Speed setting too low	— Turn speed up
Magnet jumping around erratically	— Magnet has not coupled correctly with stir bar	— Turn off and move container to recouple stirbar
Opaque fluid does not mix	— Missing stirbar	— Add missing stirbar

Disposal Responsibilities:

The equipment you purchased may contain hazardous substances that could impact the environment. Per regulations on electronic devices in the European Community, you must use the appropriate disposal systems to avoid exposure of these substances to the environment. The disposal systems will reuse or recycle hazardous materials from your equipment responsibly.

The crossed-out wheeled bin symbol invites you to use those systems.

If you need more information on the collection, reuse and recycling systems, please contact your local or regional waste administration. You can also contact us for more information on the environmental performance of our products.

Return for Repair

Important:

Transporting hazardous materials without a permit is a violation of federal law.

Heathrow Scientific LLC will not accept any return that has not been appropriately cleaned and decontaminated.

In the unlikely event of return, or when damage to the unit necessitates return, contact Heathrow Scientific LLC and obtain return authorization before sending your stirrer for service. Product must be decontaminated before transport.

Warranty

IN NO EVENT WILL HEATHROW SCIENTIFIC® LLC'S OBLIGATION UNDER THIS WARRANTY EXCEED THE PRICE OF THE PRODUCT.

Limited Warranty

Heathrow Scientific LLC warrants that Mini Stirrer will be free from defects in workmanship and materials for 5 years from the date of purchase.

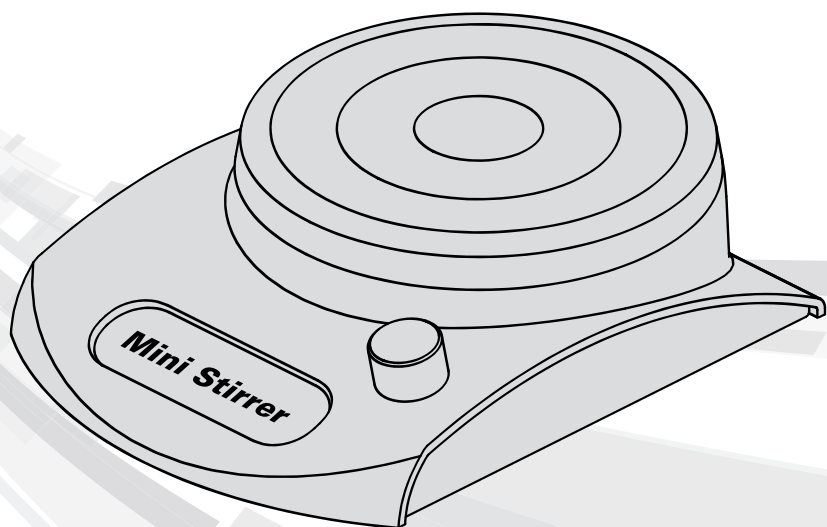
If you believe that there is a defect in the product, you must, during the warranty period, notify Heathrow Scientific LLC, provide proof of purchase, and return the product to Heathrow Scientific LLC with a Return Authorization form. To obtain a Return Authorization form, please call 1-847-816-5070. If Heathrow Scientific LLC is properly notified and, after inspection, confirms that there is a defect and the warranty period has not expired, Heathrow Scientific LLC will repair, modify, or replace the product, at its sole option, at no charge.

OTHER THAN THIS LIMITED WARRANTY, HEATHROW SCIENTIFIC LLC MAKES NO WARRANTIES, EXPRESS OR IMPLIED, WITH RESPECT TO THE QUALITY OR PERFORMANCE OF THE PRODUCT, INCLUDING BUT NOT LIMITED TO THE IMPLIED WARRANTIES OF MERCHANTABILITY AND FITNESS FOR PARTICULAR PURPOSE WHICH ARE HEREBY DISCLAIMED AND EXCLUDED. HEATHROW SCIENTIFIC LLC WILL IN NO EVENT BE LIABLE FOR ANY LOSS OF USE, LOSS OF PROFITS, CONSEQUENTIAL, SPECIAL, EXEMPLARY OR PUNITIVE DAMAGES.

THIS WARRANTY DOES NOT COVER:

- ANY DEFECT OR DAMAGE CAUSED BY IMPROPER OR UNREASONABLE USE OF THE PRODUCT. (MINI STIRRER) IS DESIGNED FOR USE ONLY BY TRAINED LABORATORY TECHNICIANS. USE BY ANYONE ELSE WILL VOID THIS WARRANTY.)
- ANY PRODUCT THAT HAS BEEN, IN HEATHROW SCIENTIFIC® LLC'S SOLE JUDGEMENT, TAMPERED WITH, ALTERED, OR REPAIRED BY ANYONE OTHER THAN HEATHROW SCIENTIFIC LLC.
- ANY PRODUCT THAT IS INOPERATIVE BECAUSE OF: (a) WEAR OCCASIONED BY USE, (b) NEGLIGENCE, (c) ACCIDENT, (d) INCORRECT MAINTENANCE, OR (e) USE UNDER ABNORMAL CONDITIONS OF TEMPERATURE, DIRT OR CORROSION, OR USE WITH ABRASIVE OR CORROSIVE MATERIALS.
- ACCESSORY PARTS, SUCH AS RUBBER PARTS, THAT ARE DAMAGED BY LIQUIDS OR MISUSE.

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All rights reserved. Printed in China.
DIE HAFTUNG UNTER DER GARANTIE VON HEATHROW SCIENTIFIC® LLC ÜBERSTEIGT UNTER KEINEN UMSTÄNDEN DEN PREIS DES PRODUKTS.
©Heathrow Scientific® LLC